




PIPETMAN P2 AND PETMAN P10

Models P2 and P10 of PIPETMAN® L, G, Neo, and Classic are equipped with a dual-position tip-ejector adapter or tip-ejector extension.

It allows the use of PIPETMAN DIAMOND® Tips D10 (short) and DL10 (long), for better precision and protection from contamination when using microtubes.

DESCRIPTION	P/N
Extension for plastic tip-ejector for PIPETMAN P2 / P10	F2070903
Dual-position adapter for stainless steel tip-ejector for PIPETMAN P2 / P10	F144879

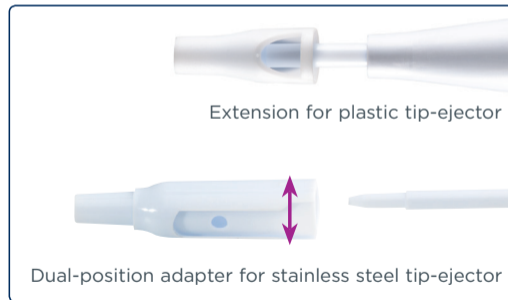


For PIPETMAN Classic, Neo, G, and L equipped with stainless steel tip-ejector

A dual-position adapter (plastic) is required to fit DL10 tips (long tips) or D10 tips (short tips).

P2 and P10 models are delivered with the adapter in place, and ready to use DL10 tips.

If D10 tips are used, the adapter must be repositioned in the shorter slot as follows:

- 1 Pull the adapter down from the stainless steel tip-ejector.
- 2 Turn the adapter 180°.
- 3 Refit the adapter so that the end of the stainless steel tip-ejector engages the shorter slot of the adapter.

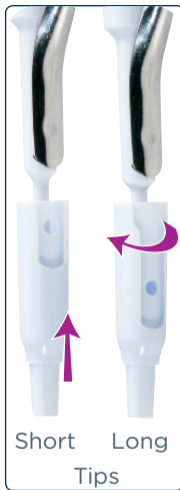

For PIPETMAN G, and L equipped with plastic tip-ejector


A tip-ejector extension is supplied to fit with D10 tips (short tips).

To fit a tip-ejector extension:

- 1 Slide the extension over the tip-holder.
- 2 Push the extension firmly onto the end of the tip-ejector until it clicks into place.

To remove a tip-ejector extension:

- 1 Gently twist the extension.
- 2 Pull it away from the pipette.

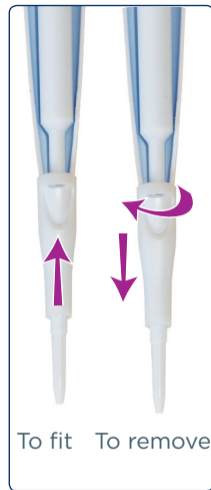

Both the dual-position adapter and tip-ejector extension are autoclavable.