



Parts

Commonly used parts (e.g., rubber inserts, TC gaskets, batteries, filters) are available from your local dealer or Drummond Scientific Company. Call us at **1-800-523-7480**.

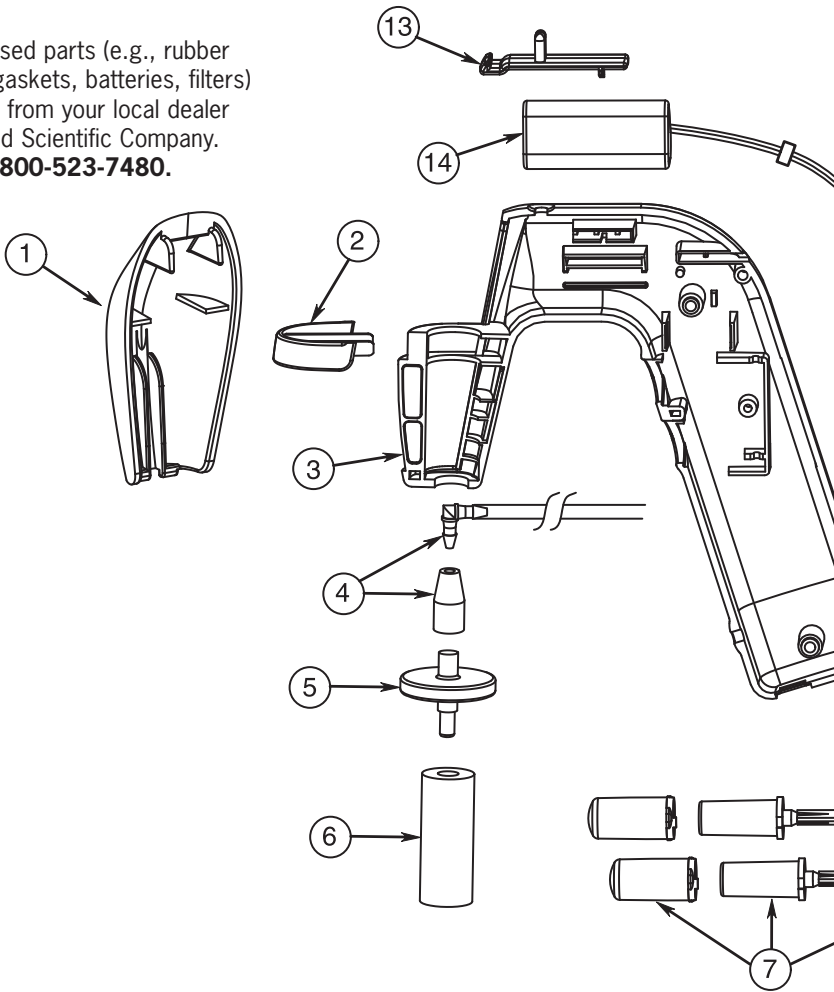

	Catalog No.	Parts Description
(1)	4-000-520	Nosepiece Cover–Front Cover–Filter Access Cover
(2)	4-000-522	Filter Lock
(3)	4-000-523	Handle Shell
(4)	4-000-524	Filter Connector–Filter/Tubing Connector
(5)	4-000-051 4-000-051-S	Self-locking Filter Self-locking Filter, Sterile (5/pk)
(6)	4-000-002	Rubber Insert
(7)	4-000-525	Button Assembly
(8)	4-000-073	XP2 Valve Body Complete w/Buttons
(9)	4-000-072	Switch Operator, XP/XL/XP2
(10)		Valve Body Slide (yellow) comes with Valve Body

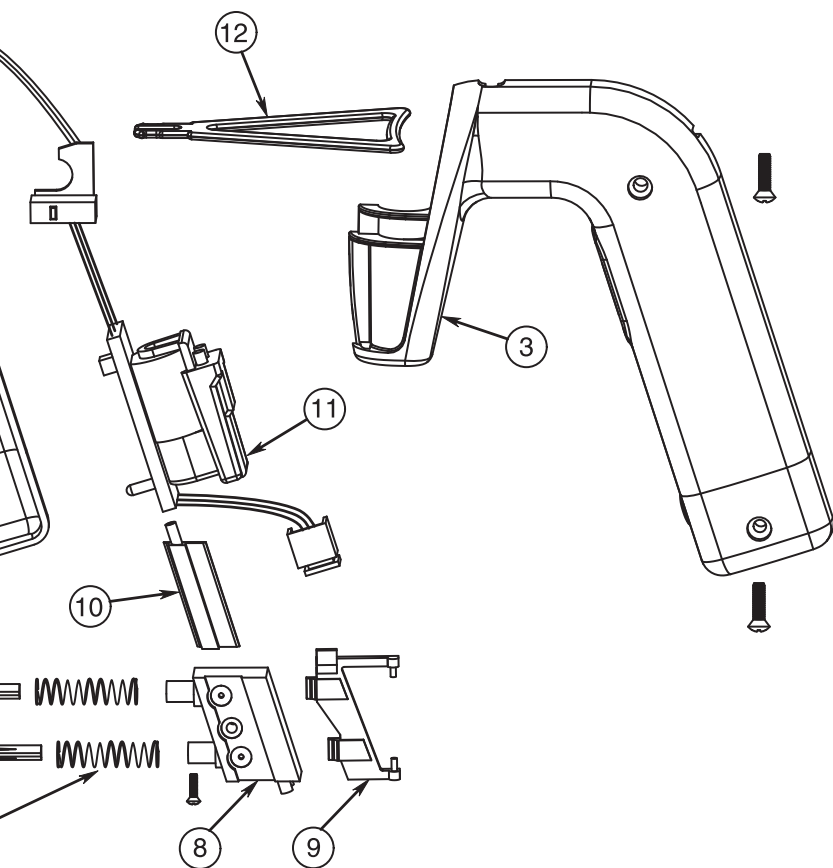

	Catalog No.	Parts Description
(11)	4-000-061	XL/XP2 Pump, Board, Wire Harness, Micro Switch with Connecting Tubing
(12)	4-000-526	XP2 Removable Stand
(13)	4-000-527	Front Cover Hold-down Clip
(14)	4-000-036	Replacement Ni-MH Battery
[Not Shown]	4-000-065	Power Supply/Charger, Input: 120V, 60Hz, 55mA Output: 6VDC, 200mA
	4-000-066	European Power Supply/Charger, Input: 230V, 50Hz, 55mA Output: DC6V, 200mA
	4-000-066-A	Australian Power Supply/Charger, Input: 240V, 50Hz, Output: 6VDC, 300mA
	4-000-066-E	UK Power Supply/Charger, Input: 240V, 50Hz, 55mA Output: DC6V, 200mA