Dispensette® HF

Mode d'emploi

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BRANDTECH
SCIENTIFIC, INC.

BrandTech Scientific, Inc.
11 Bokum Road
Essex, CT 06426-1506
USA
Tel: (860) 767-2562
Fax: (860) 767-2563
E-Mail: mail@brandtech.com
Internet: www.brandtech.com

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Safety Instructions

This instrument may be involved with hazardous materials, operations, and equipment. This Manual does not purport to address all of the safety problems associated with its use. It is the responsibility of whomever uses this instrument to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

1. **Follow the general instructions for hazard prevention and safety regulations, e.g., wear protective clothing, goggles and gloves.**
2. Observe the reagent manufacturer's information.
3. Every user must be acquainted with this Operating Manual before operation.
4. Use the instrument only for dispensing liquids, with strict regard to the defined Operating Exclusions and Limitations (⇒ pages 6-7).

- If in doubt, contact the manufacturer or supplier.
5. When dispensing, the discharge tube must always point away from the user or any other person. Avoid splashes. Only dispense into suitable vessels.
 6. Never carry the mounted instrument by the cylinder sleeve or the valve block. Breakage or loosening of the cylinder may lead to personal injury from chemicals.
 7. Never press down the piston when the screw cap is screwed on.
 8. Never remove the discharge tube or the SafetyPrime™ valve while the cylinder is filled.
 9. Never use force on the instrument.
 10. Use smooth gentle movements to operate the piston upwards and downwards.
 11. Use only original manufacturer's accessories and spare parts. Don't carry out technical modifications.
 12. Before use check the instrument for visual damages. In case of trouble (e.g., piston difficult to move, sticking valves or leakage), immediately stop dispensing. Clean the instrument according to the cleaning instructions before any further use of the instrument (⇒ pages 15-18) or contact the manufacturer.

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Application and Design

This instrument is designed for dispensing liquids, observing the following limits:

- 15 to 40 °C of instrument and reagent
- vapor pressure up to 500 mbar
- density up to 3.5 g/cm³
- viscosity up to 500 mm²/s
(dynamic viscosity [mPas] = kinematic viscosity [mm²/s] x density [g/cm³])

Materials

When the instrument is used correctly, the dispensed liquid comes into contact with only the following chemically resistant materials:

Al₂O₃, ETFE, PFA, FEP, PTFE, platinum-iridium, PP (screw cap).

Operating Exclusions

Never use this instrument for:

- liquids attacking Al₂O₃, ETFE, PFA, FEP and PTFE (e.g., dissolved sodium azide)
- liquids which are decomposed catalytically by platinum-iridium (e.g., H₂O₂)
- bases and saline solutions
- carbon disulfide
- suspensions (e.g., of charcoal) as solid particles may clog or damage the instrument.

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Operating Limitations

Liquids which form deposits may make the piston difficult to move or may cause jamming (e.g., crystallizing solutions or highly concentrated alkaline solutions).

When dispensing flammable media, make provisions to avoid static charging, e. g., do not dispense into plastic vessels; do not wipe instruments with a dry cloth.

Warning!

Never use force on a piston which is difficult to move. Immediately stop dispensing and follow the cleaning instructions (⇒ pages 15-18) or contact the manufacturer.

Unstable solutions (e.g., H₂O₂) may react with the platinum-iridium valve spring and cause changes in the reagent through catalytic action.

Compatibility of the instrument with trace material analysis must be checked by the user.

For dispensing HF:

- Observe the maximum permitted concentration of 52 %
- Always operate instrument at least once every four weeks when filled
- Only store instrument in cleaned condition (⇒ Cleaning, pages 15-18).

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Components

A	Piston seat
B	Cylinder sleeve
C	Pointer
D	Volume adjustment knob
E*	Piston
F*	Cylinder
G	Air vent cap
H*	Valve block with serial number
I*	Filling valve
J*	Filling tube
K*	Seal, PTFE
L**	SafetyPrime™ valve
M	Locking nut
N**	Backflow tube
O*	Discharge valve (integrated into discharge tube)
P*	Discharge tube
Q*	Screw cap
R*	Seal, PTFE
S*	O-Ring, Viton® (FKM), for sterile dispensing (green)
T	Mounting tool
U	Cover plate

* parts in contact with the media
 ** SafetyPrime™ valve requires backflow tube

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Attention!

The adapters supplied with the instrument are made of polypropylene (PP), and can only be applied for media which do not attack PP (⇒ page 31, ETFE-adapters).

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Warning!

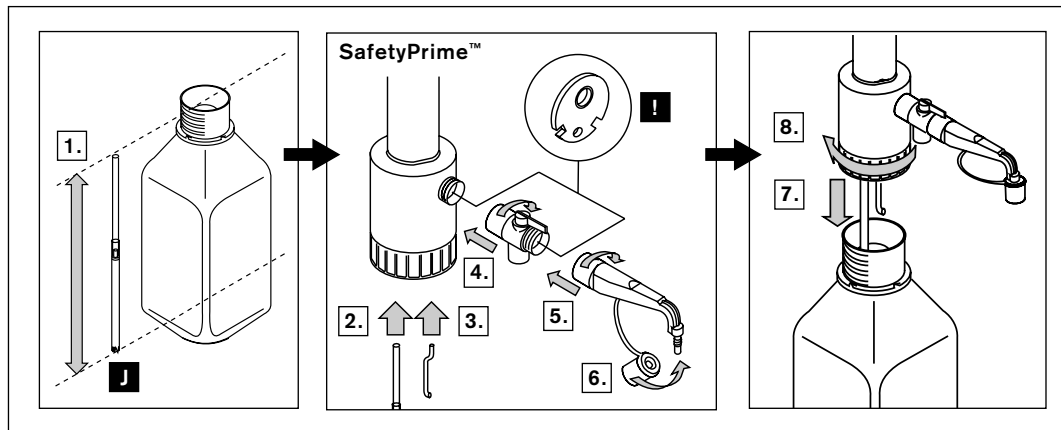
Wear protective clothing, goggles and gloves. You must follow all Safety Instructions and observe the Operating Exclusions and Limitations (⇒ pages 4-7).

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Note:

Before using the instrument for the first time, ensure it is rinsed carefully or discard first few samples dispensed.

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Attention!

To avoid splashes prime the instrument in backflow position. For small bottles use a bottle stand to prevent tipping over.

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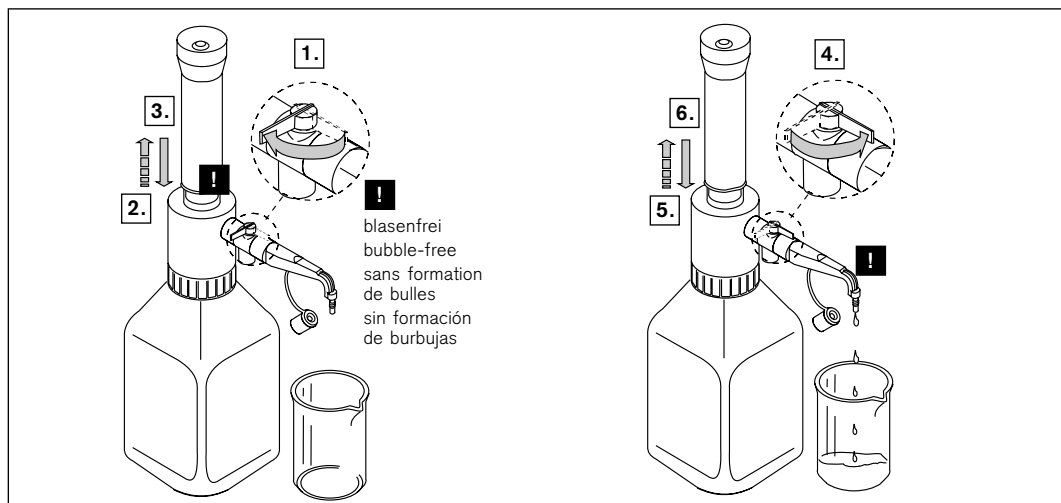
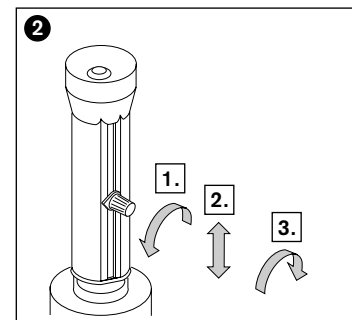
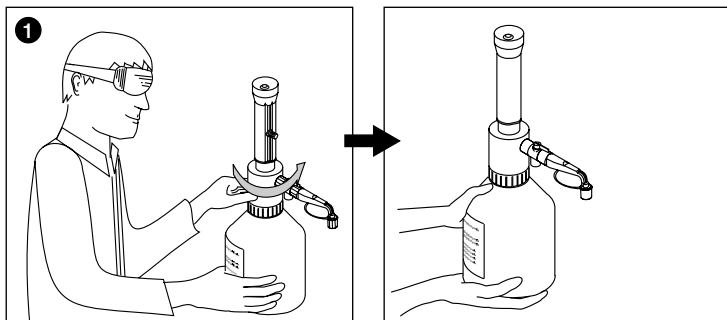
Warning!

Always wear protective gloves when touching the instrument and the bottle, especially when using dangerous liquids. Carry the mounted instrument only as shown in the figure at the right hand side.

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Warning!

Follow all Safety Instructions and observe the Operating Exclusions and Limitations (⇒ pages 4-7). The discharge tube must always point away from the user or any other person. Never press down the piston when the screw cap is pushed on. Reagent may drip off the discharge tube and screw cap.

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Attention!

After use, always leave the piston in the down position.

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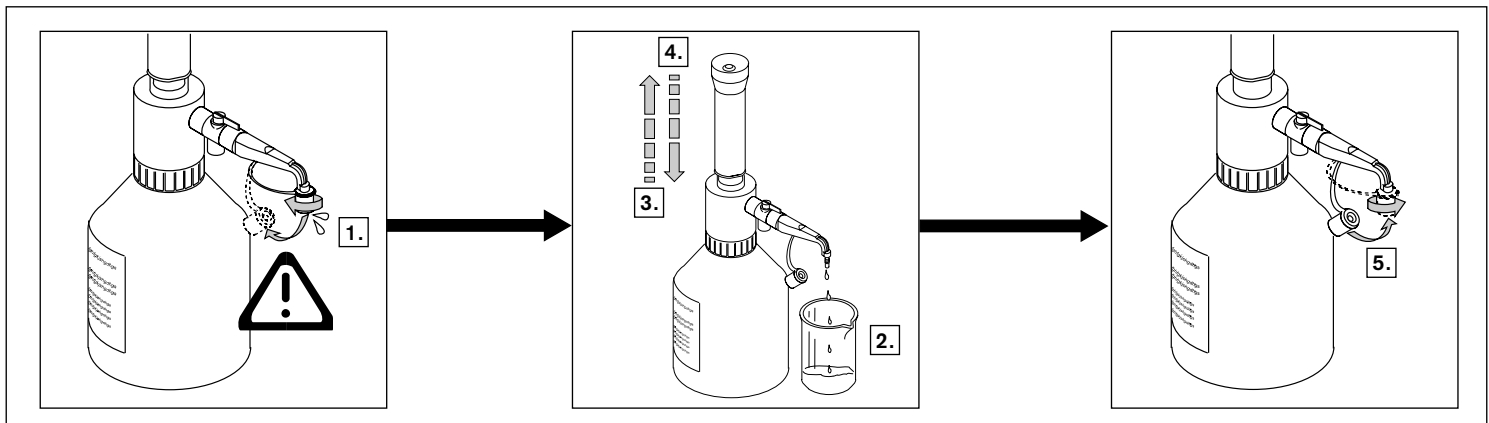
Note:

Operate instrument at least once every four weeks when filled. Only store instrument in cleaned condition (⇒ Cleaning, pages 15-18).

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Cleaning

The instrument must be cleaned in the following situations to assure correct operation:

- if the piston becomes difficult to move
- after dispensing liquids which form deposits
- before changing the reagent
- prior to long term storage
- should liquid have accumulated in the screw cap
- prior to sterilization
- before replacing the valves

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Warning!

Follow all Safety Instructions (⇒ page 4-5). The Al_2O_3 -ceramic cylinder, valves, filling and discharge tubes contain reagent. Point the valves and tube openings away from your body. Wear protective clothing, goggles and appropriate hand protection.

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Cleaning

– continued –

- ❶ Empty the instrument and unscrew.
- ❷ Mount the instrument onto another bottle filled with an appropriate cleaning solution.
- ❸ Flush the instrument.
- ❹ Empty the bottle, clean it and fill it with distilled water. Repeat step 3 and 4.
- ❺ Empty the instrument.

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
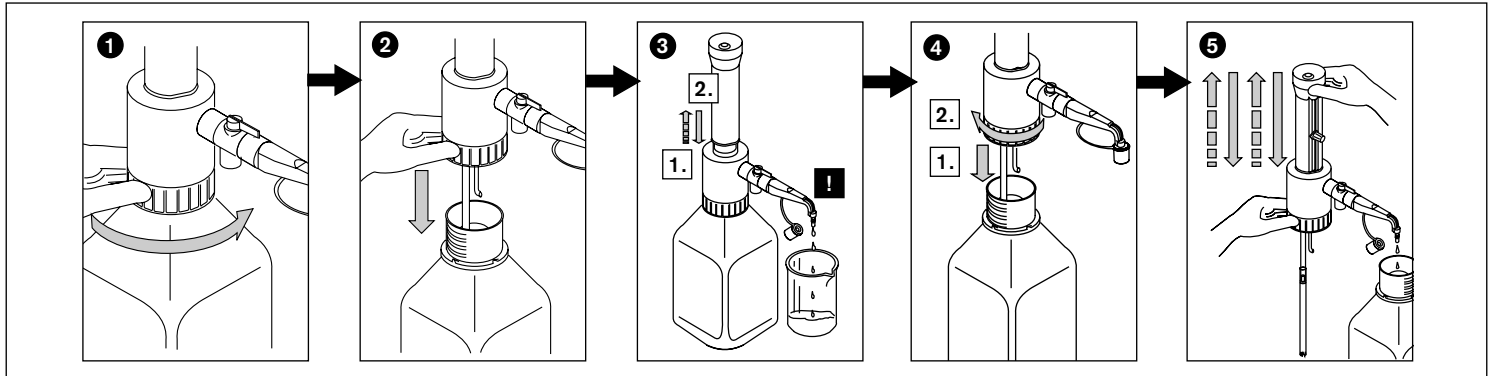
- ❻ Loosen and remove piston by pulling on the piston seat. Never exchange the piston with another instrument!
- ❼ Carefully remove deposits at the edge of the cylinder with a screw driver. Clean piston and cylinder with a bottle-brush. Flush instrument with distilled water.

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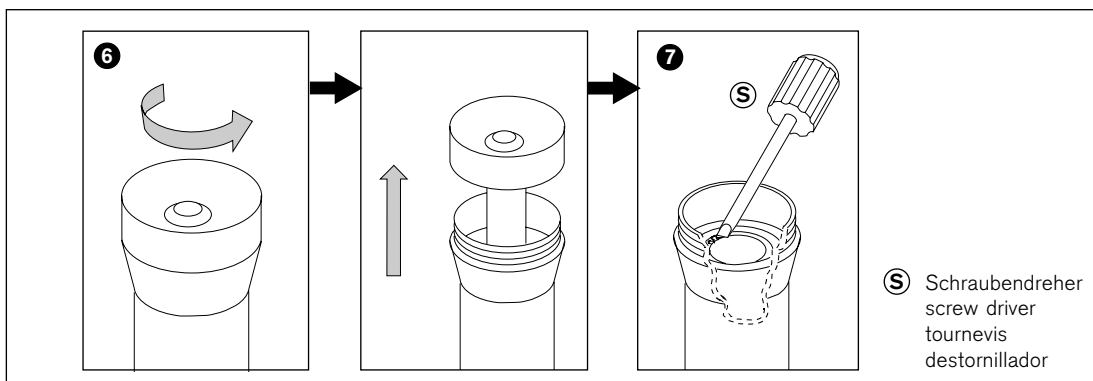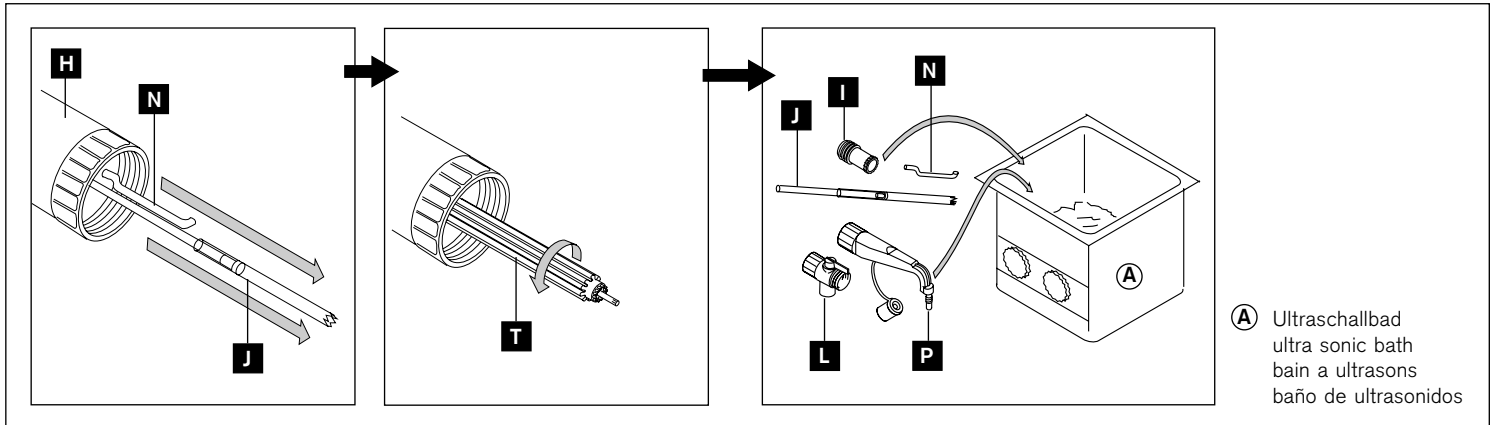
Autoclaving

This instrument withstands steam
sterilization at 121 °C (250 °F), 2 bar
absolute (15 psi) for 20 minutes in
accordance with widely used guide-
lines. It is the user's responsibility to
ensure effective autoclaving.

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**Preparation prior to the
first autoclaving**

Remove the built-in PTFE-ring (R)
and replace with the supplied elastic
O-ring (S) made of FKM.

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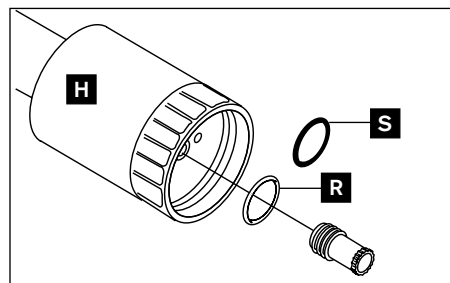
Note:

FKM has limited chemical
resistance! Do not use e.g.,
with HF!

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Autoclaving

Follow all cleaning instructions
(⇒ pages 15-18).

- 1 + 2** **Loosen** all push-in and screw connections.
- 3** **Tighten** the filling valve.
- 4** **Place** all parts (instrument with piston in down position) on a towel and autoclave. Avoid contact with metallic surfaces.

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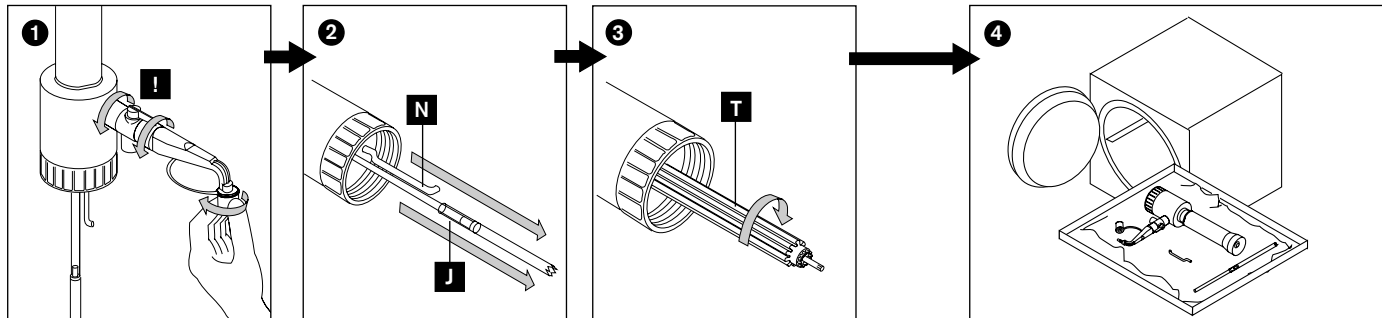
Note:

Allow to cool down in the closed autoclave to avoid deformation through too rapid cooling.
After every autoclaving, inspect all parts for deformities or damage. If necessary, replace them.
Do not reassemble the instrument until it has cooled down to room temperature.

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Checking the volume

Note:

Detailed testing instructions can be ordered from BrandTech.

- Set to the nominal volume.
- Dispense distilled water.
- Weigh the dispensed quantity on an analytical balance.
- Calculate the dispensed volume taking the temperature into account.
- Perform at least 5 dispensing and weighing operations.
- Calculate the accuracy (A%) and the coefficient of variation (CV%) by means of the formulas used in statistical quality control. The process is described, e.g., in ISO DIS 8655/2+/3. Observe the operating manual of the balance manufacturer and the corresponding standards.

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Calculations (for nominal volume)

$$\text{Mean value } \bar{x} = \frac{\sum x_i}{n}$$

x_i = results of weighings
 n = number of weighings

$$\text{Mean volume } \bar{V} = \bar{x} \cdot z$$

Z = correction factor
(e. g., 1.0029 µl/mg at 20 °C, 1013 hPa)

Accuracy

$$\text{A\%} = \frac{\bar{V} - V_0}{V_0} \cdot 100$$

V_0 = nominal volume

Standard deviation

$$s = Z \cdot \sqrt{\frac{\sum (x_i - \bar{x})^2}{n - 1}}$$

Coefficient of variation

$$\text{CV\%} = \frac{100 \cdot s}{\bar{V}}$$

Note:

For evaluation and documentation to GLP and ISO we recommend the calibrating software EASYCAL™.

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Calibration

Note:

It might be appropriate to calibrate the instrument after a longer period of usage or for specific applications.

1. Calibrate at nominal volume (⇒ page 22).
2. Calculate mean volume (result of weighing).
3. Adjust the instrument (to the calculated mean volume).

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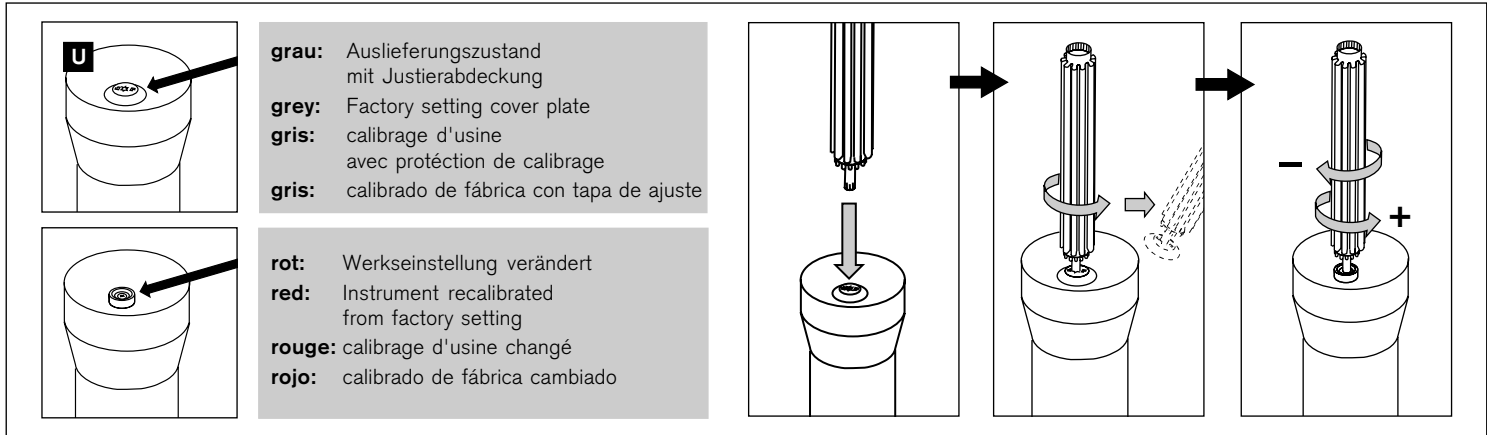
Note:

After the adjustment, it is **necessary** to make another calibration check.

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Troubleshooting

Problem	Possible cause	Corrective action
Piston is difficult to move	– Formation of crystals, dirty	– Stop dispensing immediately. Follow all cleaning instructions (⇒ pages 15-18).
Instrument does not draw in liquid	– Volume adjusted to minimum setting – Sticking filling valve	– Set to required volume (⇒ page 12). – Clean filling valve (⇒ page 18). If necessary replace the valve with sealing washer.
Instrument does not dispense liquid	– Sticking discharge valve	– Clean discharge valve.
Air bubbles in the instrument	– Reagent with high vapor pressure has been drawn in too quickly – Air not expelled from the instrument – Filling tube not firmly connected or damaged – Valves not firmly installed or damaged – Backflow tube not connected	– Slowly draw in reagent. – Remove the air (⇒ page 11). – Cleaning procedure (⇒ pages 15-18). Push the filling tube on firmly. If necessary cut off approx. 1 cm of tube at the upper end and re-connect it or replace filling tube. – Cleaning procedure (⇒ page 18). Tighten the valves using the mounting tool. If necessary, replace the valves and sealing washers. – Connect backflow tube (⇒ page 10).
Dispensed volume is too low	– Discharge tube not firmly connected or damaged – Filling tube not firmly connected or damaged – Valves not firmly connected or damaged	– Push the original manufacturer's discharge tube on firmly. Replace a deformed or damaged discharge tube. – Cleaning procedure (⇒ page 18). Push the filling tube on firmly. If this does not suffice, cut off approx. 1 cm of tube at the upper end and re-connect it or replace filling tube (⇒ page 32). – Cleaning procedure (⇒ page 18). Tighten the valves using the mounting tool. If necessary, replace valves and sealing washers.

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* Final test values (A = Accuracy, CV = Coefficient of variation) relative to the nominal capacity (maximum volume) indicated on the instrument, obtained at equal temperature (20 °C) of instrument, ambience and dist. water, and with smooth operation.

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Items supplied

1 Dispensette® HF,
1 performance certificate,
1 operating manual,
1 discharge tube,
1 filling tube (125-240 mm),
1 mounting tool,
and adapters of polypropylene of
the following size:
GL 25, GL 28, GL 32, GL 38, S 40.

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Pack of 1

Attention!

Only the following accessories
and spare parts are permitted to
use with the Dispensette® HF!

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Return for Repair

Attention:

Transporting of hazardous materials without a permit is a violation of federal law.

BrandTech Scientific, Inc. will not accept instruments that are not appropriately cleaned and decontaminated.

⇒ Therefore, contact BrandTech Scientific, Inc. and obtain return authorization **before** sending your instrument for service.

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Warranty

As provided by law, any and all warranties are null and void if the product has been misused, modified or repaired by unauthorized personnel, if the defects are caused by negligence (instruction manual, maintenance) or by normal wear and tear. Use only original manufacturer's accessory/spare parts.

For conditions and extent of warranty refer to our warranty card enclosed with the product.

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BrandTech Scientific, Inc. wird keine Geräte annehmen, die nicht ordnungsgemäß gereinigt und dekontaminiert wurden.
BrandTech Scientific, Inc. n'accepte que des appareils dûment nettoyés et décontaminés.
BrandTech Scientific, Inc. no acepta aparatos que no hayan sido debidamente limpiados y descontaminados.