GLASSMASTER™

Pipette Controller

OPERATOR INSTRUCTIONS

1 – 4

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6 – 13

RAININ
INSTRUMENT CO., INC.

Introduction

GlassMaster has been carefully designed and manufactured as a lightweight, cordless pipetting aid for use with all plastic or glass pipettes from 1–100ml.

This new instrument is constructed in recyclable materials and operates on two rechargeable, long life, environmentally friendly, 1.2V nickel metal hydride batteries. **GlassMaster** has a uniquely formed finger triggers to minimize any high pressure points. These triggers also control intake and dispense speeds.

A battery warning light (4) is triggered whenever the charge in the batteries is low. Pipetting with **GlassMaster** can also take place while the units are connected to their chargers (5).

GlassMaster has a conveniently located slide switch (3) which allows selection between three modes of operation.

HIGH Aspiration and dispense at full speed
(includes Blow-out)

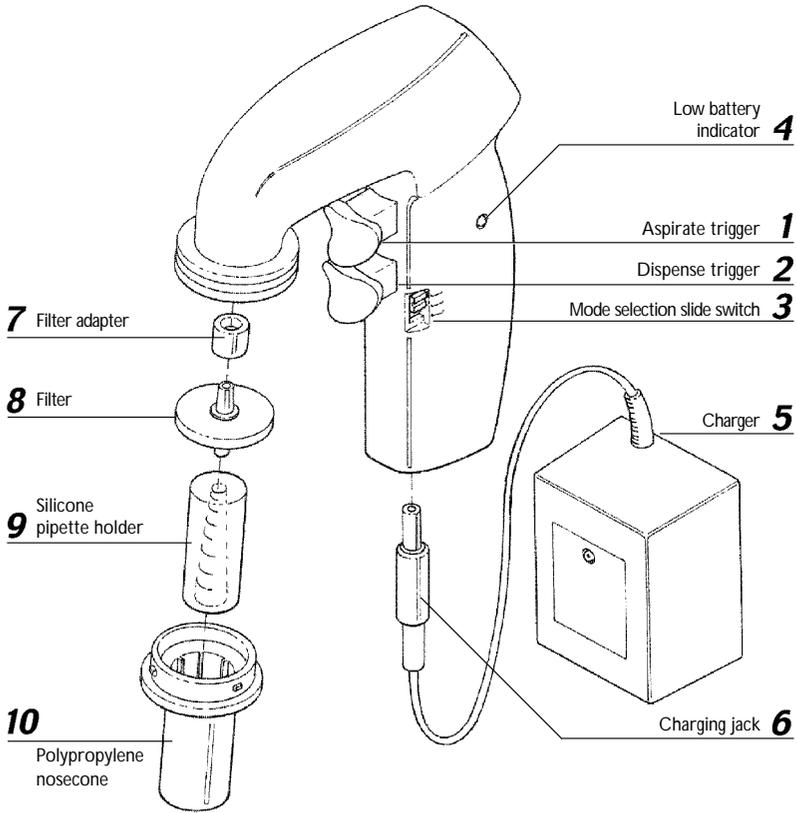
LOW Aspiration and dispense at reduced speed
(includes Blow-out)

G Aspiration at reduced speed, dispense by gravity (To Deliver)

Technical Data

Batteries	2 x 1.2V nickel metal hydride
Battery service Life	500 – 1000 cycles from empty
Charger	65mA /H for 14 hours plug-in
Overall weight	180g without stand
Filter (8)	Hydrophobic 25 mm dia 0.45 mm (fitted) plus two additional spare filters 1x 0.45 µm 1x 0.20µm
Manifold material	ABS
Case material	ABS
Nosecone material	Polypropylene (10)
Pipette holder	Silicone (9)
Pipette types	Plastic or glass 1 – 100 ml

The **GlassMaster** Pipette Controller is shipped with:

Charger; Appropriate voltage:

120V 60Hz 2-pin (USA), 230V 50Hz 2-pin (EU), 230V 50Hz 3-pin (UK)

Two spare hydrophobic filters, 1x 0.45µm and 1x 0.20µm

Bench stand, Instruction manual

Safety

GlassMaster is designed with safety in mind. Light in weight, it balances easily and safely in the hand. The triggers are carefully designed to eliminate pressure points and control of intake and delivery is therefore made easier.

When using aggressive liquids you should wear protective clothing, including goggles and gloves. If in doubt ask your safety officer. Please consult and establish appropriate health and safety practices applicable to the liquid being dispensed within your working environment.

WARNING!

- Only use this instrument within recommended limits of its chemical resistance.
- Never use **GlassMaster** with fuming liquids which are incompatible with polypropylene or silicone.
- Never use with highly flammable liquids.
- Observe the reagent manufacturer's safety information.
- Never use **GlassMaster** in an atmosphere where there is an explosion. danger.
- When not in use store on bench stand.
- Never use force when in use.
- If the instrument fails to work to specification stop pipetting. Check instructions headed Trouble Shooting (page 4).
- Only recharge with manufacturer's original charger.
- All repairs must be carried out by an authorized dealer.
- Use only the manufacturer's accessories and spare parts. See page 5.

Recharging GlassMaster

Before starting to operate carefully check mains voltage and recharge batteries.

To charge **GlassMaster**

1. Recharging temperature +10°C to +35°C
2. Plug the charger jack (6) into the base of **GlassMaster**.
3. Plug the charger (5) into a suitable wall socket. The indicator on the charger will light up.
4. Place **GlassMaster** onto bench stand
5. Allow unit to charge for 14 hours from empty to fully charged.
6. When **GlassMaster** is fully charged, the charger automatically switches to a trickle charge operation and the indicator will turn off. If the charge to **GlassMaster** is interrupted and then re-started, the charger starts its full 14 hour charging cycle again. This will not damage the battery.

Only use the charger indoors and protect against moisture.

Cleaning

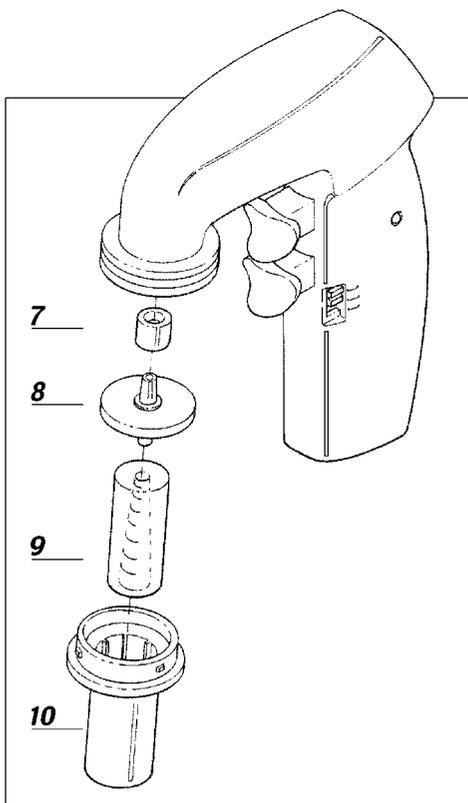
GlassMaster can be cleaned with most common laboratory disinfectants. The polypropylene nosecone (10) and silicone pipette holder (9) can be autoclaved.

To remove the nosecone and pipette holder turn the nosecone counter-clockwise. When the nosecone is removed the silicone pipette holder can also be disengaged from the filter by gently turning.

The hydrophobic filter (8) requires occasional replacement when it is wetted by the liquid and whenever suction capacity is slowing.

Replacing the Filter

The hydrophobic filter (8) will require occasional replacement, particularly when suction capacity is slowing.

Note

Only use genuine manufacturer's recommended filters and accessories as listed on page 5.

Pipetting

To aspirate

1. Carefully hold the pipette and insert it securely into the polypropylene nose cone and silicone holder.
2. Select mode of operation — High, Low, or 'G'.
3. Depress upper trigger (1) (The harder the pipetting triggers are pressed the faster the aspiration.)

To dispense

1. In BLOW-OUT mode, (mode switch high or low) depress the lower trigger (2) (The harder the pipetting triggers are pressed, the faster the dispense).
2. If 'G' gravity mode is chosen, the upper aspirate trigger automatically operates at low speed while the lower trigger opens the pipette to atmosphere, allowing liquid to run out.

WARNING!

Please ensure that no liquid enters the instrument. **GlassMaster** should never be used without the filter (8). The filter also prevents liquid from entering the instrument. Where liquid has entered please clean and replace filter.

Trouble Shooting

TROUBLE	POSSIBLE CAUSE	REMEDY
Suction capacity slow	Filter blocked, dirty or damaged	Clean or replace membrane filter
	Batteries low	Recharge
Pipette drips	Pipette not securely positioned	Re-attach in silicone pipette holder
	Pipette holder or pipette damaged	Replace

GlassMaster is virtually service-free. If the recommendations above do not resolve the problem, please contact your authorized dealer.

Note

When the meniscus level has been reached liquid should not drip out of the pipette. If the pipette does drip, check that the adaptor assembly and filter are correctly assembled and the pipette is properly located into the silicone pipette holder.

Return for Repair

In the unlikely event of repair, or where the damage to the unit necessitates return to your distributor or the manufacturer, please decontaminate the unit and complete the required decontamination certificate where appropriate. Please also give details of the fault.

Voltages

US	120V 60Hz 2-pin
Europe	230V 50Hz 2-pin
Japan	100/120V 50/60Hz 2-pin

Ordering Information

GlassMaster Pipette Controller

Cat. No.	Description
GM-100	GlassMaster Pipette Controller (1 to 100 mL)

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© 1997 Rainin Instrument Co, Inc
GlassMaster is a trademark of Rainin Instrument Co, Inc